



ImpaRATOR™

Lyophilized

STORE AT

-20°C



FOR RESEARCH USE ONLY

Instructions for Use

ImpaRATOR™ Lyophilized 2000 units (G1-IR1-020)

Process 2 mg O-glycoprotein



www.genovis.com/ifu-G1-IR1

Lyophilized Enzyme for O-glycan-specific Protein Digestion

ImpaRATOR is an O-glycan-dependent protease that catalyzes the hydrolysis of the peptide bond adjacent to glycosylated serine or threonine residues in glycoproteins and glycopeptides. It cleaves N-terminally of serine or threonine residues modified with mucin-type O-glycans, including sialylated species.

Mucin-type O-glycans are required for ImpaRATOR activity and the enzyme will not digest unmodified serine or threonine residues, or at N-glycosylation sites of glycoproteins. The enzyme accepts a broad range of O-glycan structures, including sialylated core 1 and core 2 structures as well as the Tn antigen.

ImpaRATOR is derived from *Pseudomonas aeruginosa* and expressed in *E. coli*. The enzyme contains a His-tag and has a molecular weight of 97 kDa.

UNIT DEFINITION

One unit ImpaRATOR Lyophilized digests $\geq 95\%$ of 30 μg etanercept at least at one site when incubated in TBS (50 mM Tris-HCl, 150 mM NaCl), pH 7.6 at 37°C for 30 minutes. For O-glycopeptide mapping applications, one unit ImpaRATOR Lyophilized digests 1 μg O-glycosylated protein when incubated in TBS, pH 7.6 at 37°C for 2 hours.

CONTENT AND STORAGE

ImpaRATOR Lyophilized is supplied lyophilized in TBS (50 mM Tris-HCl, 150 mM NaCl), pH 7.6, with no preservatives added. ImpaRATOR Lyophilized is shipped at ambient temperature, and should be stored at -20°C upon arrival.

After reconstitution, the ImpaRATOR enzyme is stable for at least 1 month at +4-8°C.

ImpaRATOR is for R&D use only.

QUALITY CONTROL

ImpaRATOR Lyophilized is tested to meet the specifications and lot-to-lot consistency.

ImpaRATOR Lyophilized is tested for absence of microbial contamination with blood agar plates, Sabouraud dextrose agar plates and fluid thioglycollate medium.

YOU MIGHT ALSO BE INTERESTED IN

OpeRATOR™

O-glycan specific protein digestion

OmniGLYZOR®

Hydrolysis of N- and mucin-type O-glycans

OglyZOR™

Hydrolysis of core 1 O-glycans

GlycOCATCH™

Enrichment of O-glycopeptides

O-glycan-specific Protein Digestion

PREPARATIONS

Additional Materials Required

- Digestion buffer: TBS (50 mM Tris-HCl, 150 mM NaCl), pH 7.6.¹

Sample Preparation

Prepare the glycoprotein in the digestion buffer.

The final protein concentration in the digestion reaction should be 0.1-10 mg/ml.

1. Optimization may be required if buffers other than the recommended are used. The enzyme is active at pH 5.5-8.5.

WORKFLOW

1. Prepare ImpaRATOR

1.1 Reconstitute ImpaRATOR in 50 μ l ddH₂O to 40 units/ μ l.

2. Add ImpaRATOR

2.1 Add 1 unit ImpaRATOR / 1 μ g glycoprotein.²

3. Digestion

3.1 Incubate for 2 h³ at 37°C.

2. A higher enzyme concentration may increase the digestion efficiency of individual glycoproteins. Optimization is then required.
3. Incubation times up to 18 hours may be required for complex substrates.

USA & Canada

Genovis Inc.

10919 Technology Place Suite C, San Diego, CA 92127, USA

Phone: 1-855-782-0084 (toll free)

Fax: 1-858-524-3006

EMEA & Asia

Genovis AB

Box 790, SE-220 07 Lund, Sweden

Phone: +46 46 10 12 30

Fax: +46 46 12 80 20

support@genovis.com

www.genovis.com

All rights reserved. Genovis products may be covered by one or more patents, trademarks and copyrights owned or controlled by Genovis AB. For more information about commercial rights, please contact the Genovis team at licensing@genovis.com.

Genovis products are intended for research use only. They are not intended to be used for therapeutic or diagnostic purposes in humans or animals.

All goods and services are sold subject to Genovis' General Terms and Conditions of Sale.

© Genovis AB